

For Research Use Only  
Store at Room Temperature



**INSTRUCTION**

# MANUAL

**X-Amp™ DNA Reagent**

IB47440, IB47441, IB47442

**Model Numbers:**

IB47440  
 IB47441  
 IB47442

**Quantity:**

500 µl  
 50 ml  
 100 ml

## Introduction

IBI X-Amp DNA Reagent is designed for efficient release of DNA for direct use in PCR reactions without purification. A wide variety of samples are effectively homogenized in the reagent without any pre-treatment or subsequent bind, wash or elution steps. Simply place the sample in the reagent, follow the 2-step protocol and transfer the lysate to a PCR mix.

## Advantages

- Use DNA directly in PCR reactions
- DNA purification is not required
- 15 minute 2 step protocol
- Wide variety of sample types (tissue, blood, plant, bacteria, yeast/fungus, virus)

## Applications

Direct use of DNA in PCR reactions, multiplex PCR, Real-time PCR

## Quality Control

IBI X-Amp DNA Reagent is tested on a lot-to-lot basis according to IBI's ISO-certified quality management system. DNA from a 1 mg tissue sample is lysed in X-Amp DNA Reagent. A 5 µl aliquot of lysate is added directly into a 50 µl PCR mix.

 **During operation, always wear a lab coat, disposable gloves, protective goggles or (anti-fog) procedure mask.**

## X-Amp™ DNA PCR Reagent Protocol Procedure

Sample	Procedure
Tissue	<ol style="list-style-type: none"> <li>1. Transfer 50 µl of X-Amp DNA Reagent and 1 mg of tissue to a 1.5 ml microcentrifuge tube.</li> <li>2. Incubate for 15 minutes at room temperature or 5-15 minutes at 80°C.</li> <li>3. Mix by vortex then transfer a 2-5 µl aliquot to a 20-50 µl PCR mix.</li> </ol>
Plant Tissue <sup>1</sup>	<ol style="list-style-type: none"> <li>1. Transfer 200 µl of X-Amp DNA Reagent and 5-25 mg of tissue to a 1.5 ml microcentrifuge tube.</li> <li>2. Incubate for 15 minutes at room temperature or 5-15 minutes at 80°C.</li> <li>3. Mix by vortex then transfer a 2-5 µl aliquot to a 20-50 µl PCR mix.</li> </ol>
Whole Blood, plasma, serum	<ol style="list-style-type: none"> <li>1. Transfer 100 µl of X-Amp DNA Reagent and 5-10 µl of fluid sample to a 1.5 ml microcentrifuge tube.</li> <li>2. Incubate for 15 minutes at room temperature.</li> <li>3. Mix by vortex then transfer a 2-5 µl aliquot to a 20-50 µl PCR mix.</li> </ol>
Saliva	<ol style="list-style-type: none"> <li>1. Transfer 100 µl of X-Amp DNA Reagent and 10 µl of saliva to a 1.5 ml microcentrifuge tube.</li> <li>2. Incubate for 15 minutes at room temperature or 10 minutes at 80°C.</li> <li>3. Mix by vortex then transfer a 2-5 µl aliquot to a 20-50 µl PCR mix.</li> </ol>

Bacteria <sup>2</sup>	<ol style="list-style-type: none"> <li>1. Transfer 100 µl of X-Amp DNA Reagent and 1-5 µl of bacteria culture to a 1.5 ml microcentrifuge tube.</li> <li>2. Incubate for 15 minutes at room temperature or 10-15 minutes at 80-90°C.</li> <li>3. Mix by vortex then transfer a 2-5 µl aliquot to a 20-50 µl PCR mix.</li> </ol>
-----------------------	---

NOTE: <sup>1</sup> For plant species with high levels of polysaccharide inhibitors, increase the sample amount by 2-3 times per volume of X-Amp™ DNA PCR Reagent. Plant tissue homogenization using a bead beating instrument or pestle and mortar with liquid nitrogen will facilitate DNA release.  
<sup>2</sup> E.coli can be efficiently lysed in X-Amp™ DNA PCR Reagent for 15 minutes at room temperature. However, to efficiently disrupt the bacteria cell wall of gram (+) bacteria, 3 hour incubation at room temperature or 10-15 minutes at 80°C is required.

## **IBI Stable Temp PCR Reagents – Ideal for use with IBI X-Amp DNA Reagent**

IB43101	No-Dye TAQ Master Mix	100 RXNS
IB43102	No-Dye TAQ Master Mix	500 RXNS
IB43103	No-Dye TAQ Master Mix	1000 RXNS
IB43111	TAQ HotStart No-Dye Master Mix	100 RXNS
IB43112	TAQ HotStart No-Dye Master Mix	500 RXNS
IB43113	TAQ HotStart No-Dye Master Mix	1000 RXNS
IB43120	TAQ KEEN GREEN Master Mix	100 RXNS
IB43121	TAQ KEEN GREEN Master Mix	500 RXNS
IB43122	TAQ KEEN GREEN Master Mix	1000 RXNS



SCIENTIFIC

7445 Chavenelle Road • Dubuque, IA 52002  
800-253-4942 • (563) 690-0484 • info@ibisci.com • [IBISCI.com](http://IBISCI.com)

